


Malaise Trap, Townes Style

Package Contents:

1 Malaise Trap | 1 bottle connecting ring | 1 collecting bottles

Assembly:

1. Please note that guy ropes, tent pegs or poles, support poles are not included in the trap which are otherwise require for the installation of the trap
2. It is suggested to use either wooden or aluminum poles for the support or you can find supports in the forest or field itself
3. Do not leave much gap at the bottom as the trap is made of fine quality mesh which allow easy passage of air without much resistance for proper standing of the for longer duration in the field
4. Four extra collection bottles are provided with the trap as an accessory. You may wish to carry collected insects in these bottles to a laboratory

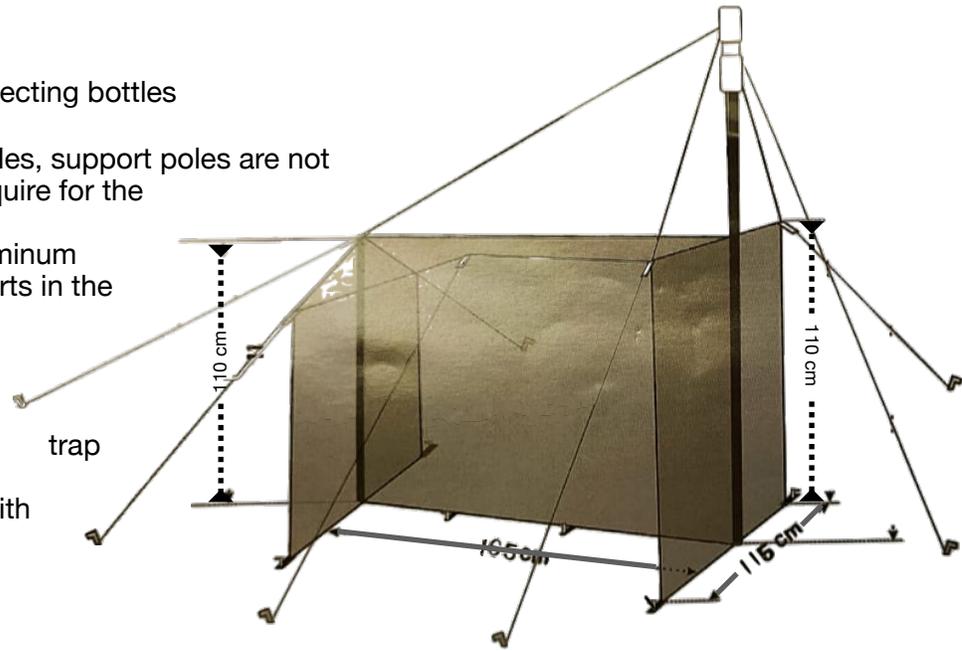

Note: Please do not over pull or push the connectors, joints or assembly poles into the provisions given in the Malaise trap. Please read instructions and understand the procedure and pitfalls before attempting to install. Wrong or improper installation attempt may damage the product. For any queries or support or suggestions please visit www.labitems.co.in or write to us at info.labitems@gmail.com


Malaise Trap, Townes Style

Package Contents:

1 Malaise Trap | 1 bottle connecting ring | 1 collecting bottles

Assembly:

1. Please note that guy ropes, tent pegs or poles, support poles are not included in the trap which are otherwise require for the installation of the trap
2. It is suggested to use either wooden or aluminum poles for the support or you can find supports in the forest or field itself
3. Do not leave much gap at the bottom as the trap is made of fine quality mesh which allow easy passage of air without much resistance for proper standing of the trap for longer duration in the field
4. Four extra collection bottles are provided with the trap as an accessory. You may wish to carry collected insects in these bottles to a laboratory

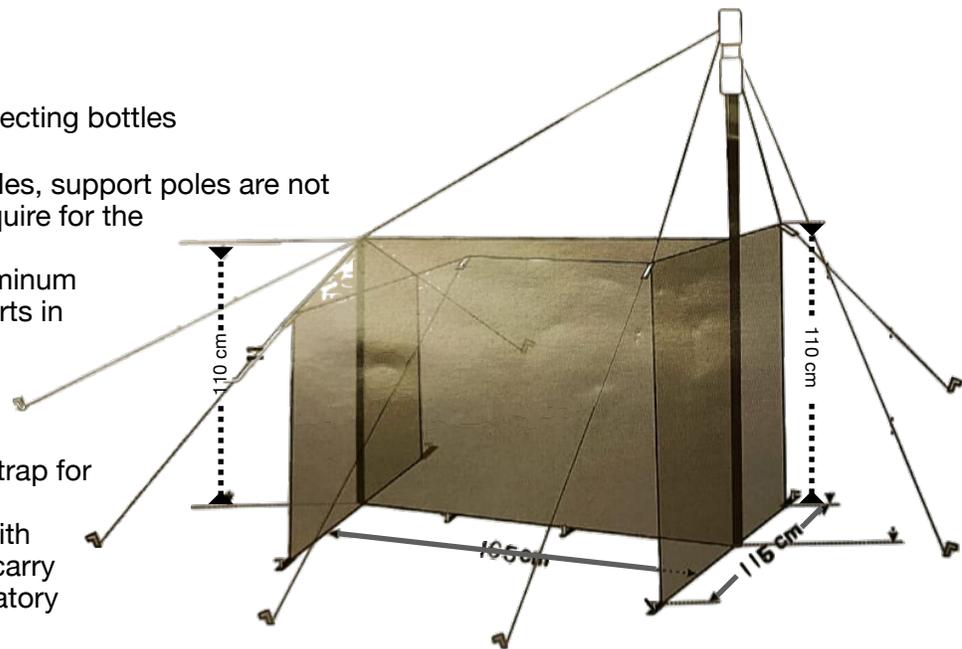

Note: Please do not over pull or push the connectors, joints or assembly poles into the provisions given in the Malaise trap. Please read instructions and understand the procedure and pitfalls before attempting to install. Wrong or improper installation attempt may damage the product. For any queries or support or suggestions please visit www.labitems.co.in or write to us at info.labitems@gmail.com